让流入人体里的制药用水最安全
MAKE PHARMACEUTICAL WATER SAFEST TO FLOW INTO THE HUMAN BODY

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Shanghai EDA Medical & Pharmaceutical Technology Co., Ltd. was established in 2011, has a rich experience team familiar with pharmaceutical industry rules. EDA can provide the EPC services to pharmaceutical fluid process system: Design, Factory modularization fabrication, Site installation, Site commissioning, Validation, Concentration automation one-station service to help the customer reduce the interface and improve the control and execution efficiency of the project execution.

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We provide integrated engineering solutions for worldwide pharmaceutical industries in compliance with EU GMP, FDA, WHO, PIC/S cGMP principle, etc. Our company cooperates closely with well-known scientific research institutions and pharmaceuticals industry designing institutes to provide reasonable project design, high-quality equipment, and customized service for different customers all over the world.

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EDA consistently holds the quality policy of "unremitting quality pursuit, everything is done for customers". As a famous machinery manufacturer, we believe that the sustainable development of our enterprise lies in the endless pursuit of perfect quality. Provided with modern production and inspection facilities, our company carries out production in strict accordance with the ISO9001 quality management system. By keeping to the operation principle of "seeking quality and pursuing perfection", EDA personnel will provide new and old customers with first class products and perfect service!

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We provide you with products are repeatedly adjusted to ensure production efficiency, reduce downtime, increase security and enhance competitive advantage. EDA carries out strict inspection on every product. The inspector is highly responsible in each component, coupled with the most advanced testing equipment and strict scientific management, so that each product can withstand the careful selection of customers.

- ① 激光切割机
Laser cutting machine
- ② 全自动成品抛光机
Automatic finished products polishing machine
- ③ 全自动抛光机
Automatic polishing machine

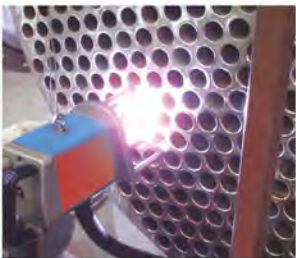
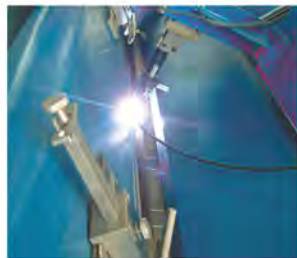
- ④ 全自动卷圆机
Automatic rounding machine
- ⑤ 全自动氩氟焊
Automatic argon fluorine welding
- ⑥ 全自动氩氟焊
Automatic argon fluorine welding

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Perfect quality relies on elaborate design and advanced equipment

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The latest digital technology is adopted for direct conversational programming and the key parts are subjected to mechanical machining such as drilling, cutting, milling and grinding to ensure the ultimate precision. We have applied brand new industrial concept and powerful technical strength to independently design and manufacture first-class products, thus completely realizing the simple and programmable operations for the consistent products.

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Strict management and elaborate manufacturing
accomplish perfect quality

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Monitoring the whole process of production is the guarantee to product quality. It is our tenet to avoid the occurrence of any defect and prevent quality problems from damaging our brand. We have established a complete set of reasonable and orderly inspection procedures. Guided by the thoughts of pursuing zero failure, We will never sell one unqualified product to clients. We believe that it not only means the responsibility for customers, but also for ourselves.

INSPECTION EQUIPMENT

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① 弹簧拉力试验
Spring tensile test

② 化学分析
Chemical analysis

③ 超声波探伤
Ultrasonic flaw detection

④ 材料光谱仪
Material spectrometer

⑤ 无损测厚仪
Nondestructive thickness gauge

⑥ 纯蒸汽品质鉴定仪
Pure steam quality identifier

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The best solution for any microbial fermentation process. Whether for single fermenters or multi-stage fermentation systems, sterile design and hygienically designed components are applied to meet the customer's highest requirements and comply with national (GMP) and international (cGMP/FDA) pharmaceutical standards.

Excellent process result in terms of quality, accuracy and repeatability which is granted by controlling pH, pO₂, temperature, agitator speed and weight through precise instrumentation.

The complete Factory Acceptance Test (FAT) in the plant enables a quick start-up at site.

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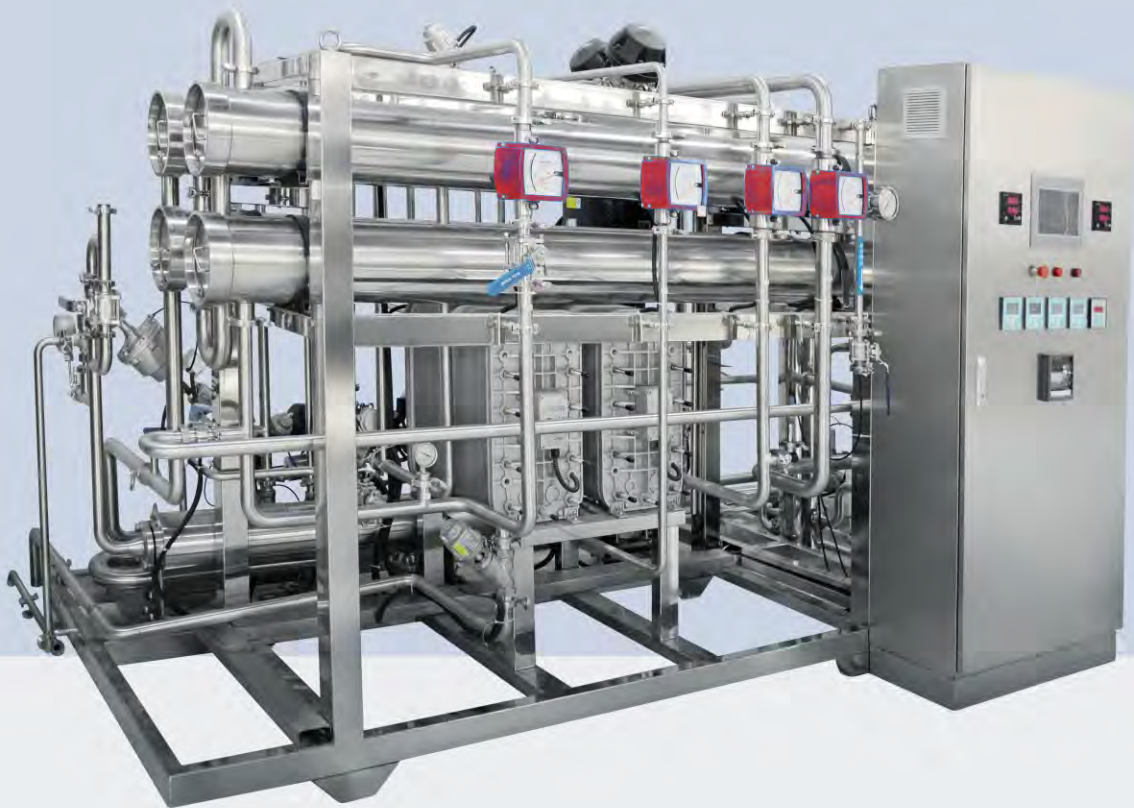
The bioreactor has the operation requirements under normal temperature and pressure to prevent bacteria and steam sterilization, and the production material is non-toxic to microorganisms or enzymes; Enzyme often because the substrate concentration is not too high, so the reactor volume is quite large; Before the reaction, it needs to access the bacteria. In order to prevent the contamination of miscellany bacteria and the decline of activity, batch reactor is generally adopted. In the fermentation process, gas, liquid and solid often coexist, and the rheological properties are complicated, so it is difficult to solve the problem of scaling up the design of bioreactor by using the principle and method of chemical reaction engineering.

Therefore, the bioreactor has the characteristics of resistance to steam sterilization, reaction is often gas, liquid, solid three-phase coexist, can prevent the contamination of bacteria and activity decline, normal temperature and pressure operation, production of tight no gap to prevent bacteria, reactor volume is quite large, non-toxic to microorganisms or enzymes of the material production and so on.

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The purified water preparation unit adopts the most advanced primary and secondary reverse osmosis + EDI process purified water treatment technology and more user-friendly 3D design, conforming to ergonomics with elegant and beautiful appearance and simple and convenient operation. It can effectively remove all kinds of salts and impurities in water, so as to make the water produced meet the quality requirements for process water in such industries as biology, medicine, medical devices, food and purified drinking water production. The system has such advantages as high automation degree, one-touch starting, stable quality of water produced, low running costs, environmental protection and convenient maintenance.

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Technical features:

1. It adopts PLC + touch screen automatic control, featured by good stability and reliability, elegant and beautiful appearance, one-touch starting, easy and convenient operation.
2. The user-friendly 3D design conforms to ergonomics. The height and position of touch screen, switch, instrument and meter are in line with average height of people, making operation easy.
3. The operations of various components such as pumps, valves and tank liquid levels are provided with dynamic monitoring pictures. (They can be upgraded into remote monitoring if necessary.)
4. It is provided with three levels of management authority such as administrator, engineer and operator.
5. It is easy to switch between manual and automatic modes. It is provided with a special alarm function to avoid misoperation by people other than operators.
6. Purified water electrical conductivity is monitored in a real-time manner. Water at the water return end is kept at a flow speed of at least 1m/s to lower too fast microbial growth at the purified water pipe network.
7. It is provided with the user-friendly function of alarm recording and indication. In case of expired replacement cycle of filter materials, water overflow, water shortage, low pressure or over pressure, such event will be recorded the touch screen event record column. In case of abnormal situations in water quality, pressure or flow, alarms will be given.
8. Each section is provided with the function of discharging nonconforming water to the previous section to ensure stable and safe water quality. (If water quality restores to be conforming in 5 minutes, the record will be made on the touch screen record column, and an alarm will be given when the time elapses.)
9. In order to realize no dead corner in all sections, if there is no fluctuation two hours after water overflow in the purified water tank, micro circulation in the whole system will be triggered to avoid microbial growth in the pipe because of no flow for a long time.
10. It is provided with the water quality correction function and the function of warning in case of minimum water quality, so that no emergency will occur.

MULTI-EFFECT WATER DISTILLATOR UNIT

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The Multi-effect water distillator unit is suitable for in vitro diagnosis, hemodialysis, injection raw water, pharmaceutical water and bioengineering laboratories in such industries as biology, medicine, medical devices and food.

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Technical features:

1. All parts that contact raw water, pure steam and water for injection are made of 316L.
2. The stainless steel surfaces are subject to electrochemical polishing to ensure the quality of injection water and improve equipment service life.
3. The external preheater, condenser and first effect evaporator adopt double tube sheet design, meeting cGMP requirements and effectively eliminating the possibility of low level media contaminating the system.
4. All hot surfaces are provided with heat insulation layers and surrounded by 304 stainless steel housings externally.
5. There is no moving part on the equipment, so its operation is quiet, safe and reliable.
6. It adopts reliable PLC control and standardized procedures.
7. It is provided with complete document verification and service system.
8. It is easy to operate and maintain.
9. It adopts PLC automatic control technology to meet different control requirements of customers. Full considerations have been given to the system, from the most basic pure steam quality judgment and distribution to higher class controls such as continuous adjustment of all media, linkage with the storage system, hierarchical password control as well as control of evaporation changes.
10. Machine performance plant test: The machine test is carried out according to the stipulated procedures in the machine test workshop by simulating the actual work environment.

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The Pure Steam Gernerator is a kind of equipment that produces pure steam by adopting deionized water as raw water and using steam heating. With such features as reasonable structure, simple operation, high thermal efficiency and low energy consumption, it is a piece of ideal equipment for sterilization in such industries as hospitals, medicine, electronics, food and beverage, and a preferred device that meets cGMP requirements in the pharmaceutical industry.

The Pure Steam Gernerator consists of evaporator, preheater and electrical automatic control part. In order to ensure pure steam quality, the evaporator adopts a unique three-level separation method: that is, falling film evaporation, gravity settling and screen defoaming.

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The automatic dispensing unit is used for culture medium preparation for cell culture and microbial culture, and online automatic preparation and CIP/SIP for the dispensing system of injection powder and liquid injection production lines. According to process requirements, it is equipped with operating platform, pipe, valve, instrument and meter, electric part and control software, which form complete modular equipment for completing FAT, SAT and other services. The equipment is manufactured in strict accordance with GMP and GAMP5, and DQ, IQ, OQ protocols are established according to user requirements. Tests are carried out in strict accordance with FAT, SAT requirements, and we can help users perform PQ.

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CIP/SIP system is one of the prerequisites of food Pharmaceutical production sanitary standards. It can remove the active mixed pollution and insoluble particles of impurity, reduce or remove the micro chemicals and heat source that pollute the productions. GMP recommends that food and pharmaceutical companies use the CIP online cleaning system and SIP online sterilization system, which limits related industries because of the high expense in overseas completely automatically CIP/SIP system. The specialists in our company import and digest international advanced technique, and successfully make automatical CIP/SIP system with high automation, simple operation, and reasonable price.

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System composition

The intelligent CIP system is a centralized system, which consists of injection water tank, CIP tank, double tube sheet heat exchanger, liquid delivery pump, return pump, temperature controller, liquid level indicator, electrical conductivity detector, PLC touch screen, control cabinet and system valve and pipe fittings.

Biopharmacy: used for CIP/SIP in equipment and pipes for culture medium preparation, fermentation, harvest, ultrafiltration, purification, inactivation and material transfer

Aseptic preparations: used for CIP/SIP in equipment and pipes for material preparation, material transfer and sterile connection

Aseptic powders and injection: used for CIP/SIP in equipment and pipes for material preparation, material transfer and sterile connection

Sterile bulk drugs: used for CIP/SIP in equipment and pipes for material preparation, material transfer and sterile connection
It improves process traceability, effective validates and revalidates reliability of production equipment and reduces production risks.

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According to URS, we provide overall material distribution solutions, clean pipes, valves, and instruments and meters. The sanitary dead corner, CIP/SIP, electric control and site construction quality can meet requirements for new GMP and EU GMP certification. In addition, we supply full sets of qualification documents.

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LIST OF CUSTOMERS

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衷心感谢广大合作伙伴对上海俄达的支持与厚爱，未来，上海俄达将竭诚为广大用户提供更优质的产品，更周全的服务。一起共谋发展，共图未来！
We cordially thank our partners who have always given support and favor to EDA . In the future, EDA will sincerely provide customers with higher quality products and more considerate services. Let's develop together for a resplendent future.

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